

visicolor® ECO Hydrazin

Hydrazine / Idrazina

436 nm


Method(e) / Método

5301


0.05–0.40 mg/L N₂H₄

Test 5-30


REF 931 230

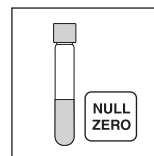


Probe / sample
échantillon / muestra

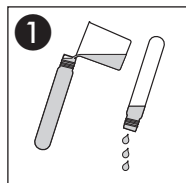


Küvetten spülen
Rinse test tubes
Rincer les éprouvettes
Lavar los tubos

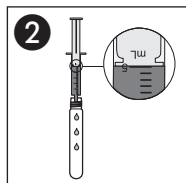


5 mL Probe
Sample
Echantillon
Muestra

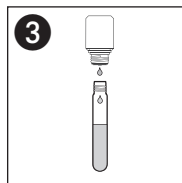


5 ∆ N₂H₄-1

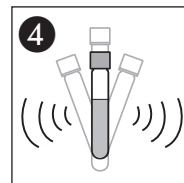


Schütteln
Shake
Agiter
Agitar

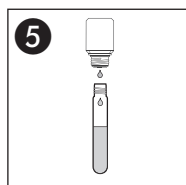


5 ∆ N₂H₄-2

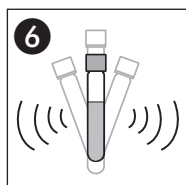


Schütteln
Shake
Agiter
Agitar

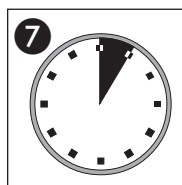

5'00 min

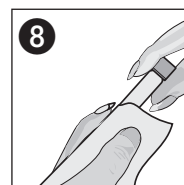

Säubern
Clean
Nettoyer
Limpiar

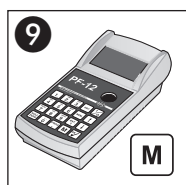

Messung
Measurement
Mesure
Medición



Meerwasser
Sea water
Eau de mer
Agua de mar