

# visicolor® ECO Sulfid

## Sulphide / Sulfure / Sulfuro

620 nm


Method(e) / Método

5941


0.05–0.80 mg/L S<sup>2-</sup>

### Test 5-94


REF 931 294

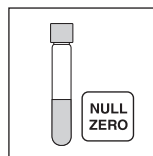


Probe / sample  
échantillon / muestra

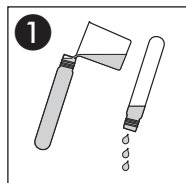


1  
Küvetten spülen  
Rinse test tubes  
Rincer les éprouvettes  
Lavar los tubos

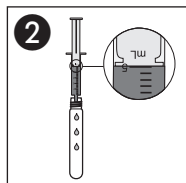


2  
5 mL Probe  
Sample  
Echantillon  
Muestra

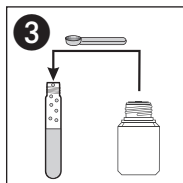


3  
1 mL S-1

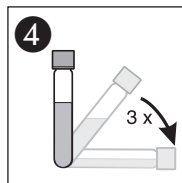


4  
Umschwenken  
Shake gently  
Secouer légèrement  
Mezclar volteándolo

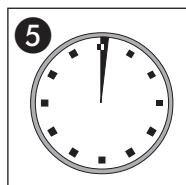


5  
1'00 min

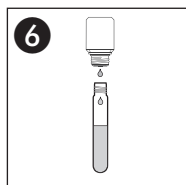


6  
5 mL S-2

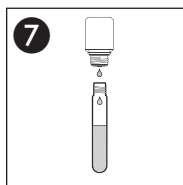


7  
5 mL S-3

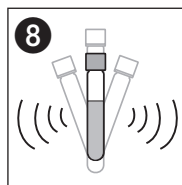


8  
Schütteln  
Shake  
Agiter  
Agitar

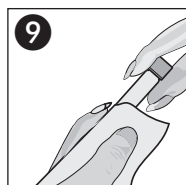

9  
Säubern  
Clean  
Nettoyer  
Limpiar

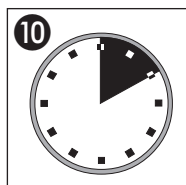

10  
10'00 min

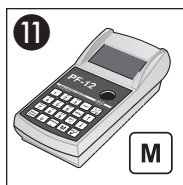

11  
Messung  
Measurement  
Mesure  
Medición



Meerwasser  
Sea water  
Eau de mer  
Agua de mar