




10  
Thermokappe aufschrauben  
Close with thermo cap  
Fermer avec thermo cap  
Cerrar con thermo cap

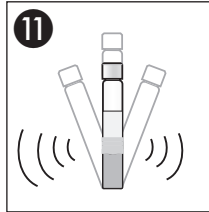


11  
Schütteln  
Shake  
Agiter  
Agitar

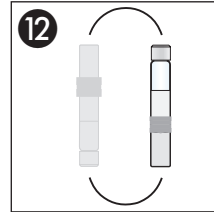


12  
Küvette um 180° drehen  
Turn tube upside down  
Retourner tube  
Girar tubo

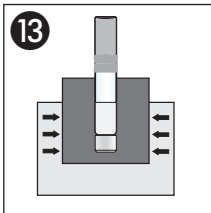


13  
120 °C / 2 h

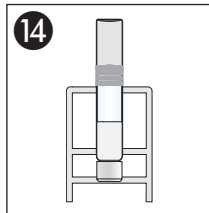


14  
60 min abkühlen lassen  
allow to cool  
laisser refroidir  
dejar enfriar

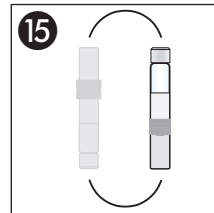

15  
Küvette um 180° drehen  
Turn tube upside down  
Retourner tube  
Girar tubo

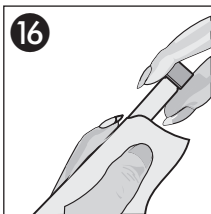

16  
Säuber  
Clean  
Nettoyer  
Limpiar

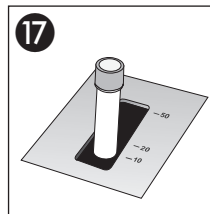

17  
Messung  
Measurement  
Mesure  
Medición